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E-mail: bioprep1@otenet.gr www.bioprepare.gr

**TECHNICAL DATA SHEET****PRODUCT: CHROMagar™ Mastitis****REFERENCE: 020366**Date 1st Edition:

7th 2009

Date 4th Revision:

6th 2024

DESCRIPTION

CHROMagar™ Mastitis is a selective chromogenic medium in a bipartite (split) Petri dish for the isolation and differentiation of pathogenic Gram(+) cocci and Gram(-) bacteria that cause mastitis.

PRINCIPLE OF THE METHOD

CHROMagar Mastitis is a new tool for the rapid and simple differentiation of the bacteria causing the infection. This new product consists of two different media: one for Gram(+) bacteria and the other for Gram(-). CHROMagar™ Mastitis GP isolates and differentiates Gram(+) cocci. CHROMagar™ Mastitis GN isolates and differentiates Gram(-) bacteria.

FORMULA	g/litre
CHROMagar™ Mastitis GP	
Chromogenic and selective mix	4,4
Peptone, yeast and extract	20,0
Salts	5,0
Agar	15,0
GP supplement	2ml

Appearance: Light beige, opaque

Final pH 6.9 ± 0.2 at 25 °C.

FORMULA	g/litre
CHROMagar™ Mastitis GN	
Chromogenic and selective mix	1,2
Peptone, yeast and extract	17,0
Agar	15,0

Appearance: Light beige, opaque

Final pH 7.0 ± 0.2 at 25 °C.**PRECAUTIONS**

CHROMagar™ Mastitis is an in-vitro laboratory diagnostic material and should only be handled by qualified people in the laboratory. This material contains peptones and extracts of animal origin. The certificates regarding the origin and health status of the animals do not fully guarantee the absence of transmissible pathogens. For this reason, it is recommended that these materials be treated as potentially infectious, with the usual safety precautions (avoiding ingestion or inhalation). Plates should always be handled with gloves and in Laminar flow Class II, to avoid contamination mainly by saprophytic fungi. If the plate is cracked or the bag has a hole, do not use it. Do not use petri dishes if there are signs of microbial contamination. The thickness of the agar must be 4 - 5 mm and the material without cracks, dryness or other signs of deterioration. After the expiry date the material is unfit for use. In case of contact with the skin, wash immediately with plenty of water and soap. Positive samples must be destroyed according to the hygienic rules prescribed for the management of contaminated samples.

STORAGE CONDITIONS

The Petri dishes must be stored at **2–12 °C** in their original packaging until use. Prolonged storage at temperatures below **2 °C** generates excess moisture within the medium, increasing the risk of contamination. Freezing, even momentarily, destroys the medium. Excessive heating should also be avoided. The dishes can be used until the expiration date indicated on the label. After opening the airtight packaging, any unused dishes should be stored in the same bag until their expiration date. For transportation, our stability studies have shown that the plates can remain at **6–25 °C** for up to **4 days** or at **25–40 °C** for **24 hours** without compromising product performance.

INSTRUCTIONS FOR USE**Presence/Absence Testing:**

The infectious sample should be delivered to the laboratory promptly, without delay, and protected from extreme heat or cold. If processing is delayed, the sample should be inoculated into an appropriate transport medium and kept at 2–12 °C until plated. Allow the plates to reach room temperature before use. The agar surface must be dry before inoculation. Inoculate the sample at one edge of the plate, then streak across the surface in parallel lines using a loop to isolate individual colonies. Incubate aerobically at 35–37 °C for 20 to 28 hours.

Bacterial Enumeration:

Perform a 1:10 dilution of the sample using an appropriate diluent (e.g., MRD).

Inoculate 0.1 ml onto the plate and spread over the entire surface.

Incubate at 35–37 °C for 24–48 hours under aerobic conditions.

Count the colonies and multiply by 100 to calculate the number of bacteria per 1 ml of sample.

If bacterial count is too high, perform further dilutions (1:100, 1:1000, etc.).

INTERPRETATION OF RESULTS – CHROMagar™ Mastitis GP

Streptococcus agalactiae forms light blue colonies.

Streptococcus uberis forms metallic blue colonies.

Staphylococcus aureus forms pink colonies.

Gram-negative bacteria do not grow.

INTERPRETATION OF RESULTS – CHROMagar™ Mastitis GN

Gram-negative bacteria:

Escherichia coli → red to brick-red colonies

Klebsiella, *Enterobacter*, *Citrobacter*, *Serratia* → blue colonies, some with pink halos

Proteus, *Morganella*, *Providencia* → colorless to beige colonies with brown halos

Proteus vulgaris → blue colonies with brown halos

Pseudomonas → yellow colonies with a green tint in the medium

LIMITATIONS OF THE CHROMagar™ Mastitis GP METHOD


Some *Staphylococcus aureus* strains may produce colonies that are pale pink to colorless. Therefore, confirmation with Coagulase and other biochemical tests is necessary.

LIMITATIONS OF THE CHROMagar™ Mastitis GN METHOD

Some *E. coli* strains may produce colonies that are pale pink to colorless. Final identification may require additional testing such as biochemical or immunological tests (e.g., latex agglutination).

QUALITY CONTROL

CHROMagar™ Mastitis GP	Colony Characteristics
<i>Streptococcus agalactiae</i>	Blue-green
<i>Streptococcus uberis</i>	Metallic blue
<i>Staphylococcus aureus</i>	Pink with halo
Gram(-) Βακτήρια	Inhibited
CHROMagar™ Mastitis GN	
<i>Escherichia coli</i>	Pink to red
<i>Klebsiella</i> , <i>Enterobacter</i> , <i>Citrobacter</i> .	Metallic blue
<i>Proteus mirabilis</i>	Brown with halo
<i>Pseudomonas</i>	Cream, translucent
<i>Candida albicans</i>	White, opaque, small colonies
Gram (+) κόκκοι	Inhibited

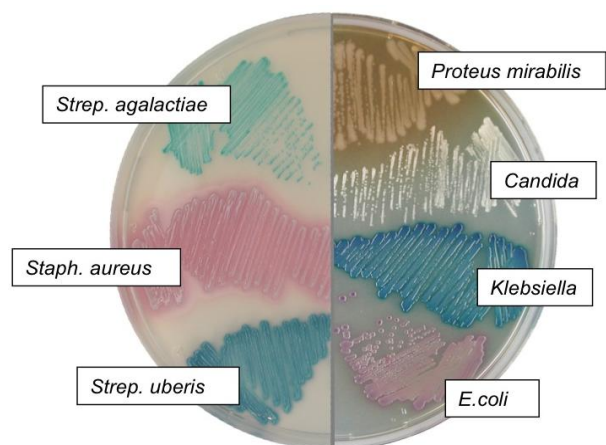WASTE DISPOSAL OF WASTE

Materials that show no growth can be considered as non-hazardous waste and disposed of accordingly. Materials that show colony growth must be disposed of according to the guidelines for infectious or potentially infectious waste. The laboratory is responsible for the proper management of infectious waste according to its nature and level of risk and must handle and dispose of it (or assign its management and disposal) in compliance with the applicable regulations.

SPECIFICATIONS

CHROMagar™ MASTITIS - CE

ITEM	CODE	PACKAGE	STORAGE	SHELF LIFE
Split Petri Dish 9 cm	020366	10 pieces	2 – 12 °C	3 months

Produced in Greece by the company Bioprep, in accordance with Regulation (EU) 2017/746. Basic UDI-DI: 5212037714010402WH. EDMA Code: (14 01 04 02) Chromogenic ID Media (Plates). Bioprep is certified according to the standards: EN ISO 9001:2015 / ELOT EN ISO 13485:2016 / DY8d/1348/2004.

REFERENCES

(1) O. Gaillot et al. 2000. Evaluation of CHROMagar Staph aureus, a New Chromogenic Medium, for Isolation and Presumptive Identification of Staphylococcus aureus from Human Clinical Specimens. JCM 38, N°4: 1587-1591. (2) A. Carricajo et al 2001. Performance of the Chromogenic Medium CHROMagar Staph aureus and the Staphychrom Coagulase Test in the Detection and Identification of Staphylococcus aureus in Clinical Samples J.C.M. 39: 2581-2583. (3) Z. Samra et al. 2004. Optimal detection of Staphylococcus aureus from clinical

IN VITRO MANUFACTURER'S DATA

G. PAPANIKOLAOU & CO

PRODUCTION LABORATORIES OF CULTURE MEDIA

Potamou 5, Industrial Area Keratea, Attica

P.O. Box: 4893, Postal Code: 9001 - Tel: +30 2299066113. Fax: +30 2299066112

E-mail: bioprep1@otenet.gr www.bioprep.gr