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CHROMagar Candida, a new differential isolation medium for the presumptive identification of clinically important *Candida* species.

Odds F. et al.

1994, Journal of Clinical Microbiology, 32 : 1923-1929. Use of CHROMagar Candida medium for isolation of yeasts from dental samples.

Beighton D. et al.

1995. Journal of Clinical Microbiology, 33 : 3025-3027.

Application of CHROMagar Candida for rapid screening of clinical specimens for *Candida albicans*, *Candida tropicalis*, *Candida krusei*, and *Candida (Torulopsis) glabrata*.

Pfaller M.A. et al.

1996. Journal of Clinical Microbiology, 34 : 58-61. Routine use of CHROMagar Candida medium for presumptive identification of *Candida* yeast species and detection of mixed fungal populations.

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E-mail: bioprep1@otenet.gr www.bioprepare.gr

DESCRIPTION

The Contact Plate Chromagar Candida is a specialized contact plate used for the isolation, cultivation, and identification of fungi of the Candida genus, and it is recommended for microbiological surface monitoring.

PRINCIPLE OF THE METHOD

In the contact plate (contact plate), the agar material can come into direct contact with the test surface. With the Contact Plate Chromagar Candida, the user can monitor surfaces such as benches, walls, product conveyor belts in industry, production machinery, tools, etc. This method using contact plates is recommended by ISO standards 14698-1 and 18593. Meat peptones provide nitrogen, vitamins, minerals, and amino acids essential for growth. Yeast extract is a source of vitamins, particularly B-group vitamins. Sodium chloride provides essential electrolytes for transport and regulates the osmotic balance of the medium. Chloramphenicol is a broad-spectrum antibiotic that inhibits a wide range of Gram-positive and Gram-negative bacteria. Bacteriological agar is the solidifying agent.

COMPOSITION	g/litre
Peptone	10.2
Chromogenic mix	22.0
Chloramphenicol	0.5
Agar	15,0

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STORAGE AND TRANSPORT CONDITIONS

The petri dishes must be stored at **2–12 °C** in their packaging until the time of use. Prolonged storage at temperatures below **2 °C** creates excess moisture inside the material, posing a risk of contamination. Freezing, even momentarily, destroys the material. Excessive heating should also be avoided. The petri dishes can be used until the expiration date indicated on the label. For transport, our stability studies have shown that the dishes can remain at **18–25 °C** for **10 days** or at **25–40 °C** for **48 hours** without affecting the product's performance.

INSTRUCTIONS FOR USE

Place the entire surface of the medium onto the surface you wish to test, which must be completely dry. Press lightly for 8–12 seconds. Place the petri dish in the incubator with the lid facing down. Incubate as indicated in the table below.

SURFACE	INCUBATION TIME	TEMPERATURE (With increased humidity)
Environmental surfaces	1–4 weeks	25 +/- 2 °C
Human monitoring	1–4 weeks	25 - 30 °C

RESULTS

After incubation, count the colonies assuming the surface area of the petri dish is 25 cm². The levels of microorganisms depend on the type of surface. Below is a table with an example for food surfaces:

Microorganisms	Interpretation
One colony / 25 cm ²	Excellent
2–10 colonies / 25 cm ²	Good
11 or more colonies / 25 cm ²	The surface must be cleaned

INTERPRETATION OF RESULTS

Candida albicans forms green-colored colonies.

Candida tropicalis forms colonies with a metallic blue color.

Candida krusei forms fuzzy pink colonies.

Candida kefyr & *Candida glabrata* form colonies ranging from purple to brown.

Other *Candida* strains form colonies with colors ranging from white to pink.


LIMITATIONS OF THE METHOD

Some bacteria may be quite fastidious and may grow very poorly or not at all.

GENERAL CHARACTERISTICS OF QUALITY CONTROL


Microorganism	Colony Growth / Color
<i>Candida albicans</i> ATCC® 10231	Green (2–6 mm – Incubation 24–48 hours)
<i>Candida tropicalis</i> NCPF3111	Metallic blue (1–2 mm – Incubation 48–72 hours)
<i>Candida krusei</i> NCPF3100	Fuzzy pink (4–10 mm – Incubation 24–48 hours)
<i>Candida glabrata</i> NCPF3112	Purple to brown (2–3 mm – Incubation 24–48 hours)

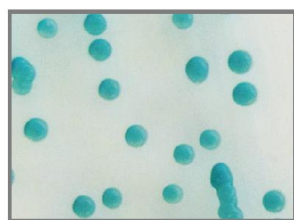
C.albicans

C.krusei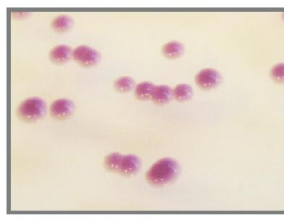

C.glabrata

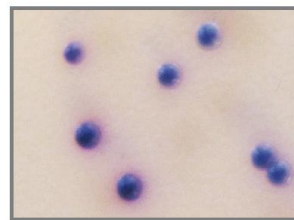C.tropicalis

DISPOSAL OF MATERIALS IN WASTE

Materials that show no growth can be considered non-hazardous waste and disposed of accordingly. Materials that show colony growth must be disposed of in accordance with guidelines for infectious or potentially infectious waste. The laboratory is responsible for the proper management of infectious waste according to its nature and level of risk and must manage and dispose of (or assign the management and disposal of) it according to the applicable regulations.

SPECIFICATIONS

CONTACT PLATE CHROMagar™ CANDIDA - CE

ITEM	CODE	PACKAGE	STORAGE	SHELF LIFE
Petri dish 65mm	20028	10 pieces	2 – 12 °C	3 months

Produced in Greece by the company Bioprep, in accordance with Regulation (EU) 2017/746. BASIC UDI-DI: 5212037714030202WV EDMA: (14 03 02 02) Chromogenic Media for Yeasts and Fungi. The company Bioprep is certified according to the standards: EN ISO 9001:2015 / ELOT EN ISO 13485:2016 / DY8d/1348/2004.

REFERENCES

- CHROMagar Candida, a new differential isolation medium for the presumptive identification of clinically important *Candida* species.
Odds F. et al.
1994. Journal of Clinical Microbiology, 32 : 1923-1929. Use of CHROMagar Candida medium for isolation of yeasts from dental samples.
Beighton D. et al.
1995. Journal of Clinical Microbiology, 33 : 3025-3027.
Application of CHROMagar Candida for rapid screening of clinical specimens for *Candida albicans*, *Candida tropicalis*, *Candida krusei*, and *Candida (Torulopsis) glabrata*.
Pfaller M.A. et al.
1996. Journal of Clinical Microbiology, 34 : 58-61. Routine use of CHROMagar Candida medium for presumptive identification of *Candida* yeast species and detection of mixed fungal populations.
Bouchara J.P. et al.
1996. Technical Report, 202-208.

IN VITRO MANUFACTURER'S DATA

G. PAPANIKOLAOU & CO

PRODUCTION LABORATORIES OF CULTURE MEDIA

Potamou 5, Industrial Area Keratea, Attica

P.O. Box: 4893, Postal Code: 9001 - Tel: +30 2299066113. Fax: +30 2299066112

E-mail: bioprep1@otenet.gr

www.bioprep.gr